R1	1 × 30 mL - Buffer pH7,6 / NADP 77mg / ATP 185mg
R2	1 × 0,6 mL - 6-PGDH 105U
R3	1 × 0,6 mL - GK 16U
C	1 × 2 mL - Control solution / Solution de contrôle

### Sample preparation:

The concentration of D-Gluconic acid in the sample used in the assay procedure has to be between 0.05 to 1.0 g/l.

### Precision:

Under the described conditions, measurement accuracy is 5% on a control solution

### Assay procedure:

Wavelength: 340nm / Optical path: 1cm / Temperature: 20-37°  
Measurement: against water or air.

	Blank	Sample
R 1	1,0 ml	1,0 ml
Water	2,0 ml	1,9 ml
Sample	0	0,1 ml
R 2	0,02 ml	0,02 ml
Mix and read DO	<b>DO 1</b>	<b>DO 1</b>
R 3	0,02 ml	0,02 ml
Mix and read Do after 15 min	<b>DO 2</b>	<b>DO 2</b>

### Calculation

Determine the values for blank and samples:

$$\Delta DO = [ DO 2 - DO 1 ]_{\text{sample}} - [ DO 2 - DO 1 ]_{\text{blank}}$$

D-Gluconic acid concentration is given by:

$$C = \frac{V \times MW}{\epsilon \times l \times v \times 1000} \times \Delta DO \quad (\text{g/L})$$

In the assay procedure:

$$C = 0,9530 \times \Delta DO \quad [\text{g/l of D-Gluconic acid in the sample}]$$

Dilution factor of the sample has to be considered in the calculation.

### Storage instructions and reagent stability

The reagents are stable up to the indicated month of expiry, if stored at 2-8 °C, contamination is avoided.

### Warnings and precautions

Do not swallow. Avoid contact with the skin and mucous membranes. Take necessary precautions for the use of laboratory reagents.

### General information on sample preparation

- Use liquid samples, clear, colorless and with pH practically 7.6, directly, or after dilution.
- Crush or homogenize solid or semi-solid samples
- Deproteinize samples containing proteins with perchloric acid; alternatively clarify with Carrez reagents.
- Extract samples containing fat with hot water

### Assay control

D-Gluconic acid reagents must be validated with the use of the standard included in the kit. The standard is ready-to-use.

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