R1	1 × 30 mL - Tampon pH9 / Buffer pH9
R2	1 × 6 mL - NAD
R3	1 × 0,65 mL - ALDH (acetaldehyde dehydrogenase)
C	1 × 0,5 g - Ethanal

Sample preparation:

The concentration of Ethanal in the sample used in the assay procedure has to be between 0.010 to 0.200 g/l.
Use clear, colorless and practically neutral liquid samples directly, or after dilution. Filter or centrifuge turbid solutions.
Ethanal is extremely volatile (boiling point = 21°C) :
- It is necessary to close tightly all containers with ethanal
- It is necessary to pipette solutions under the surface of the ethanal containing solutions

Precision:

Under the described conditions, measurement accuracy is 5% on a control solution

Assay procedure:

Wavelength: 340nm / Optical path: 1cm / Temperature: 20-37°C
Measurement: against water or air.

	Blank	Sample
R 1	1,0 ml	1,0 ml
R 2	0,2 ml	0,2 ml
Water	1,0 ml	0,9 ml
Sample	0	0,1 ml
Mix and read the DO	DO 1	DO 1
R 3	0,02 ml	0,02 ml
Mix and read DO after 10 min	DO 2	DO 2

Calculation

Determine the values for samples:

$$\Delta DO = [DO 2 - DO 1]_{\text{sample}} - [DO 2 - DO 1]_{\text{blank}}$$

Ethanal concentration is given by:

$$C = \frac{V \times MW}{\epsilon \times l \times v \times 1000} \times \Delta DO \text{ Ethanal (g/L)}$$

In the assay procedure:

$$C = 0,15722 \times \Delta DO \text{ [g/l of Ethanal in the sample]}$$

Dilution factor of the sample has to be considered in the calculation.

Storage instructions and reagent stability

The reagents are stable up to the indicated month of expiry, if stored at 2-8 °C, contamination is avoided.

Warnings and precautions

Do not swallow. Avoid contact with the skin and mucous membranes.
Take necessary precautions for the use of laboratory reagents.

Assay control

Ethanal reagents must be validated with the use of the standard included in the kit.

Prepare one control at 0,200 g/L of ethanal.

Weight 32,0 mg of ethanal, and dissolve it in 100 ml water.

Solution is stable one day. Tightly close the flask.

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