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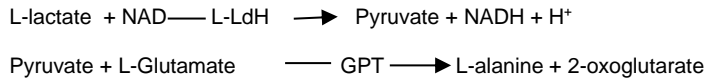
RA	1 × 100 mL – Glycyl-glycine pH10 / Glutamic acid / PVP
RB	1 × 20 mL – NAD / Glutamate Pyruvate Transaminase / L-lactate dehydrogenase / Stabilizers

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The reagent is ready to use.

Stability of reagents: see the expire date of the kit.

**Automates use:**

**Range of measurement:** The test was developed to determine the concentration in L-lactic acid ranging between 0.1 and 3 g/L in the sample with the use of the automate parameters setting supplied by Biosentec.

**Procedure of test:** For a use on automat, applications are available on request.

Follow the analytical procedure:

R1 = 250 / Sample = 3  
Incubation 2 min – Lecture 1  
R2 = 50  
Incubation 5 min – Lecture 2  
Multipoint calibration – Linear  
End-point

Wavelength: I = 340 nm / II = 700 nm

Depending on the used matrix, the calibration could be adjusted.

**Manual use:**

Contact us to get assay protocol.

**Storage instructions and reagent stability**

The reagents are stable up to the indicated month of expiry, if stored at 2-8 °C, contamination is avoided.

Onboard stability: 5 days

**Warnings and precautions**

Do not swallow. Avoid contact with the skin and mucous membranes. See the material security data sheet for more information.

**Assay control**

L-lactic acid reagents must be validated with the use of a control, available on catalogue.

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