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Cat. No. 210

210-23031

The Acids/Sugars multi solution is a titrated solution for Acetic acid, Malic acid and Lactic acid, and Glucose and Fructose. This Multi solution can be used as a Calibrator during the calibration of sequential automatic analyzer or can be used to carry out quality control.

Composition

The Multi solution is a synthetic solution and contains components suitable for carry out calibration or quality control on automatic analyzer.

Values

The reference value and measurement incertitude of the concentration of the reference solution is given on the certificate of analyse and on the label of the bottle.

Target Values	
Acetic acid	1.00 g/L
Malic acid	4.00 g/L
Lactic acid	3.00 g/L
Glucose-Fructose	8.00 g/L

The determinations were made following a validated internal method using the principle of the OIV reference methods.

Preparation and Use

Solution is ready to use.
Use the Multi solution as a calibrator or as a quality control solution.

Storage instructions and stability

The solution is stable up to the indicated month of expiry, if stored at 2-8 °C.
Onboard stability: 1 day. Recommendation is to change solution every day on the analyzer.

Warnings and precautions

Do not swallow.
Avoid contact with the skin and mucous membranes.
See the material security data sheet for more information.

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